

## How to Clean Up a Formaldehyde Spill



### Step 1: Supplies

- Mask
- Goggles
- Gloves
- Broom
- Plastic Dustpan
- Amphomag

While you may be able to clean the spill without protection if none is available, it again may be better to call in a team to avoid exposure depending on your level of training.



### Step 2: Clean Up

1. Ventilate the area by opening any windows and doors or using a fan.
2. Put on protective clothing and gear if available before getting near the spill.
3. Pour an ample amount of Amphomag onto the formaldehyde.
4. Continue to slowly add Amphomag until it completely absorbs the spill, reducing the harmful vapors and smell in the area.



### Step 3: Disposal

The formaldehyde spill is now ready to be disposed of in accordance to your local and state regulations. Use a broom and plastic dust pan to sweep up the Amphomag and dispose.