


## How to Clean Up a Hydrochloric Acid Spill

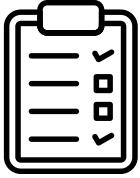

### Step 1: Supplies

Here are the supplies you will need to clean up a hydrochloric acid spill:

- Mask
- Goggles
- Gloves
- Broom
- Plastic Dustpan
- Amphomag

While you may be able to clean the spill without protection if none is available, it again may be better to call in a team to avoid acid exposure depending on your level of training.



### Step 2: Clean Up

Follow these 3 easy steps to cleaning up a hydrochloric acid spill using Amphomag.

1. Pour Amphomag slowly and directly onto the hydrochloric acid spill to contain and neutralize.
2. The Amphomag will initially turn reddish/pink to indicate that it is an acid and when fully neutralized will turn to a yellow/green color.
3. If there is any hydrochloric acid remaining anywhere, repeat step 2 until all hydrochloric acid is gone.



### Step 3: Disposal

The contained and neutralized spill is now ready to be disposed of as a non-hazardous material. Use a broom and plastic dust pan to sweep up the Amphomag.