


How to Clean Up Hydrofluoric Acid Spill

Step 1: Supplies

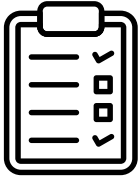

Here are the supplies you will need to clean up a hydrofluoric acid spill:

- Mask
- Goggles
- Gloves
- Broom
- Plastic Dustpan
- Amphomag

While you may be able to clean the spill without protection if none is available, it again may be better to call in a team to avoid acid exposure depending on your level of training.



Step 2: Clean Up

1. Pour Amphomag slowly and directly onto the spill to contain and neutralize.
2. The Amphomag will initially turn reddish/pink to indicate that it is an acid and when fully neutralized will turn to a yellow/green color.
3. If there is any hydrofluoric acid remaining anywhere, repeat step 2 until all hydrofluoric acid is gone.



Step 3: Disposal

The contained and neutralized hydrofluoric spill is now ready to be disposed of as a non-hazardous material. Use a broom and plastic dust pan to sweep up the Amphomag and dispose in landfill.