

## How to Clean Up Sulfuric Acid Spill



### Step 1: Supplies

- Mask
- Goggles
- Gloves
- Acid-Resistant Apron
- Broom
- Dustpan
- Water
- Amphomag



### Step 2: Clean Up

1. Check yourself to see if any of the acid got on you – or anyone else who is around – and wash off any acid found.
2. Open the windows and doors and wear protective clothing if available.
3. Pour an ample amount of Amphomag on the sulfuric acid. Amphomaga will turn red to indicate an acid.
4. Amphomag will turn a yellow/green color when it has been fully neutralized.



### Step 3: Disposal

Once Amphomag has neutralized the sulfuric acid spill, it can be swept up and disposed.